


UHPLC Analysis of Antifungals on Titan™ C18

The extremely rapid (~3 minutes) resolution of six antifungal compounds on an Titan C18 column is shown in this application. Titan particles feature totally porous, monodisperse silica. Fluka LC-MS Ultra CHROMASOLV solvents were used to supply low background interference and low particulate contaminants for robust, trouble-free operation. Cerilliant CRMs provided reliable identification and quantification.

market focus	Pharmaceutical (small molecule)
column	Titan C18, 5 cm x 3.0 mm I.D., 1.9 µm particles (577126-U)
mobile phase	[A] 5 mM ammonium acetate pH 6.8 unadjusted in 80:20, water:acetonitrile; [B] 5 mM ammonium acetate pH 6.8 unadjusted in 5:95, water:acetonitrile;
gradient	0% B; to 100% B in 3 min; 100% B held for 0.6 min
flow rate	1.0 mL/min
pressure	4350 psi (300 bar)
column temp.	35 °C
detector	UV, 205 nm
injection	1 µL
sample	6 antifungal 50 µg/mL in 70:30, water:methanol
Application No.	G006255

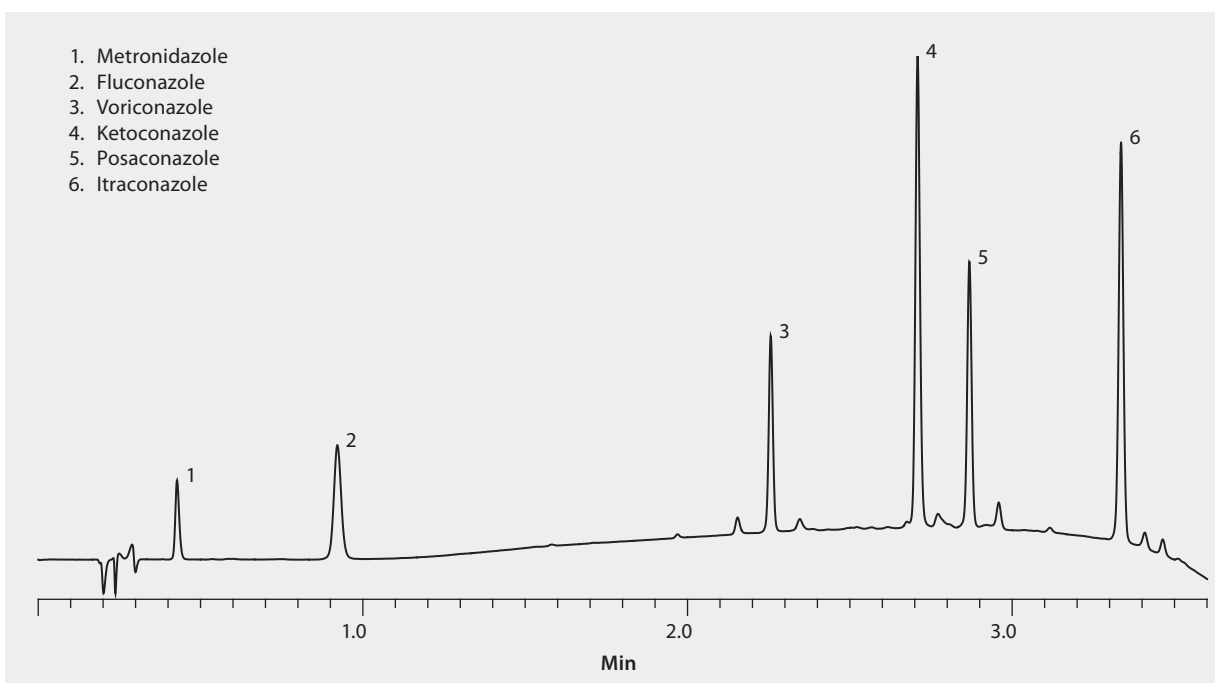

Related Products

analytical column

Titan™ C18 UHPLC Column, 1.9 micron (Supelco 577126-U)

mobile phase component

Acetonitrile (Fluka 14261)

Ammonium acetate (Fluka 14267)

Water (Fluka 14263)

standard

Fluconazole solution (Cerilliant F-031)

Itraconazole solution (Cerilliant I-015)

Ketoconazole solution (Cerilliant K-004)

Metronidazole solution (Cerilliant M-183)

Posaconazole solution (Cerilliant P-103)

Voriconazole solution (Cerilliant V-032)